User Guide NUL-205

English

Español

NeuLog logger sensors and modules

NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

Note: NeuLog products are intended for educational use.

NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at www.NeuLog.com. Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

Note: There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to www.NeuLog.com before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

Quick start procedure

PC or Mac Computer – Wired Connection

Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at www.NeuLog.com as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

Procedure:

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

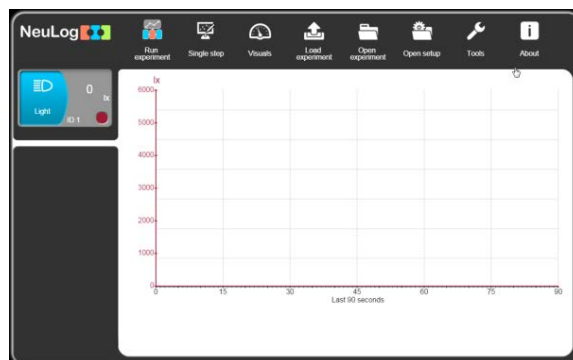6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- 'Run Experiment' with the connected sensors.
- 'Single Step' experiment.
- 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- 'Load Experiment' from the sensors' flash memory.
- 'Open Experiment' from a saved file.
- 'Tools' for changing the software settings or sensor ID.

Tablet, smart phone or any device with WIFI connectivity

Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

Procedure:

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: www.NeuLog.com. Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
 - a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

Tablet, smart phone with Bluetooth connectivity

Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

Procedure:

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

Operation with NeuLog Graphic Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

Procedure:

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

Off-line experiments

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

Materials needed:

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

Materials needed to configure your offline experiment:

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

Procedure:

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

Operation with NeuLog Digital Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

Procedure:

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

Note: If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

NeuLog Oxygen Logger Sensor Guide NUL-205

NeuLog oxygen logger sensor NUL-205

The NeuLog oxygen sensor can be used for any science experiment where oxygen levels, dissolved or gaseous, are required. It is used in the fields of Chemistry, Physiology, Exercise Science, Biochemistry, Biology, Earth Science, Environmental Science, etc.

The sensor comes pre-calibrated so you can start experimentation right out of the box using this guide.

Among hundreds of possible experiments that can be performed with the NUL-205 sensor are: monitoring of ecological systems, combustion reactions, and oxygen consumption rates, studying photosynthesis, testing chemical solutions and hydrogen peroxide decomposition, and oxygen solubility.

This sensor is practical for testing in school laboratories as well as in the field and has two modes of operation: open air mode to test gaseous oxygen concentration (percentage of oxygen in the air), and dissolved oxygen mode for the testing of liquids (mg/L and percentage of dissolved oxygen in a liquid).

The measurement units for this sensor are:

- Oxygen % in air: The concentration of gaseous oxygen present in the sample.
- mg/L: Total mass of dissolved oxygen per liter of a given liquid sample.
- Oxygen % in water: The percentage of dissolved oxygen present versus the maximum level of oxygen saturation.

Sensor's usage and offset

Prior to use:

1. Remove the blue rubber cap from the NeuLog oxygen probe.
2. Unscrew the tip-end membrane cap of the probe, inside you will see a metal electrode.
3. Add 1 mL (half of the cap) of the "DO Filling Solution" directly into the tip-end which was just unscrewed.
4. Carefully reattach the tip-end of the probe by screwing it back into the probe (do not over tighten).
5. Do not replace the blue rubber cap until after testing.

Although the NeuLog oxygen sensor does not need to be calibrated prior to each use, it is recommended to offset its probe.

Open air mode:

Offset of the probe assumes a standard atmospheric oxygen concentration of 20.9% (assuming stable levels in the Earth's atmosphere at sea level).

1. Connect the oxygen sensor to a voltage source (USB-200, BAT-200, WIFI-201).
2. Wait about 5 minutes after connecting the sensor to the voltage source before you offset the sensor. For a more accurate measurement, look at the module window and wait for the reading to stabilize (it could take a little longer).
3. Press and hold the oxygen sensor's push button for 3 seconds.

Dissolved oxygen mode:

Inserting air into water with an air pump or a 50 ml syringe will saturate the water to 100% of its capacity in the specific temperature of water you are working in (using an air pump will get more accurate readings).

1. Connect the oxygen sensor to a voltage source (USB-200, BAT-200, WIFI-201).
2. Submerge the sensor in water. Make sure that the internal metal thermistor that is 4.5 cm above the oxygen probe's tip is also submerged in the water. Use water with the same temperature for offsetting and for further readings (oxygen is less soluble in high temperatures).
3. Wait about 5 minutes after connecting the sensor to the voltage source before you offset the sensor. For a more accurate measurement, look at the module window and wait for the reading to stabilize (it could take a little longer).
4. Saturate the water with oxygen using an air pump or a 50 ml syringe.
5. If you are in the "mg/L" mode, wait until you do not see any change in the oxygen concentration. This means that you pumped the maximum amount of air into the water (in the specific temperature you are working in).

6. Press and hold the oxygen sensor's push button for 3 seconds. You will see that the oxygen concentration that appears on the module window changes to the concentration that correlates to the water temperature.
7. If you are in the "% in dissolved" mode, wait until you do not see any change in the oxygen percentage. This means that you pumped the maximum amount of air into the water (in the specific temperature you are working in).
8. Press and hold the oxygen sensor's push button for 3 seconds. You will see that the oxygen percentage that appears on the module window changes to 100%.
9. When the oxygen concentration/percentage has stabilized, it means that all excess oxygen has come out of the water (as shown in the following graph).

Included with sensor

- NeuLog General Guide.
- NeuLog oxygen sensor probe attached to the sensor's body by a durable rubber-coated wire.
- 30 mL oxygen sensor filling solution (1 M KCl).
- Replacement membrane cap for the oxygen sensor.

Sensor specifications			
	In air	Dissolved (%)	Dissolved (mg/L)
Range and operation modes	0 to 25%	0 to 125%	0 to 12.5 mg/L
ADC resolution	13 bit		
Resolution	0.1%	0.1%	0.01 mg/L
Max sample rate (S/sec)	100		

Experiment Duration: 1 second to 31 days.

Sensor's features

- Fully digital data.
- Rugged plastic ergonomic case.
- Pushbutton switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).
- Attached probe cased in durable plastic.
- Thermistor for temperature compensation.

Note: NeuLog products are intended for educational use.

Videos and experiment examples

- Videos, literature and other probes can be found at www.neulog.com.
- In order to access the oxygen sensor's page, choose "Products" on the main menu and then "Oxygen logger sensor".
- In order to access the oxygen sensor's experiments, choose "Example Labs":

Respiration (B-3)

Combustion (C-5)

Gas Solubility (C-6)


Technical background

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (for example: °C, °F, Lux, %, ppm).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update. In spite of that, the probe needs to be offset prior to initial use. Please review the "Sensor offset" guide at the beginning of this document for detailed instructions.

The sensor is based on a Clark type probe that includes a platinum electrode and an AgCl electrode placed in a KCl electrolyte; all is covered by a Teflon membrane.

Oxygen molecules diffuse through the membrane oxidizing the AgCl electrode; this accepts positive potential while the electrolyte accepts negative potential according to the following equations:

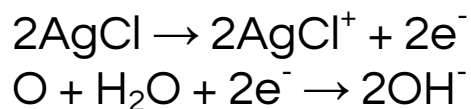The potential difference of the electrodes is converted into an oxygen measurement by the sensor's controller using the temperature compensation calculation.

The sensor offset process also depends on the internal temperature readings.

Maintenance and storage

Short-term storage (less than 24 hours):

For short term storage, fill a beaker with enough distilled water to cover the membrane cap of the oxygen probe. With the membrane cap still attached to the probe, place its tip side down in the distilled water. This easy method is ideal for short term storage, between trials for up to 24 hours.

Long-term storage (more than 24 hours):

1. Unscrew the membrane cap from the oxygen probe.
2. Discard the DO filling solution from the membrane cap.
3. Rinse the cap with distilled water and gently shake to air dry.
4. Rinse and dry the exposed metal electrode with distilled water; it can be dried with lab wipes.
5. For storage purposes: Loosely reattach the membrane cap to the oxygen probe.

Maintenance:

- Never submerge the NeuLog sensor's plastic body in any liquid.
- Do not allow liquid into the NeuLog plastic body.
- After using the probe, wipe off all excess material, liquid or residue from the oxygen probe body. Do not leave the membrane cap with filling solution for more than two days.
- Store in a box at room temperature out of direct sunlight.
- Do not put the plastic case or cord of the oxygen probe in an acid or base.
- Store the oxygen sensor's probe upright with the blue rubber tip pointing downward.
- Do not use the sensor in acid or alkaline solutions with a molarity greater than 1.0.

Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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