User Guide

NUL-209

English

Español

NeuLog logger sensors and modules

NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

Note: NeuLog products are intended for educational use.

NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at www.NeuLog.com. Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

Note: There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to www.NeuLog.com before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

Quick start procedure

PC or Mac Computer – Wired Connection

Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at www.NeuLog.com as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

Procedure:

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

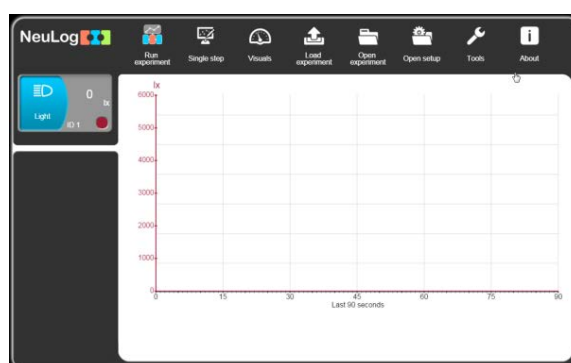6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- a. 'Run Experiment' with the connected sensors.
- b. 'Single Step' experiment.
- c. 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- d. 'Load Experiment' from the sensors' flash memory.
- e. 'Open Experiment' from a saved file.
- f. 'Tools' for changing the software settings or sensor ID.

Tablet, smart phone or any device with WIFI connectivity

Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

Procedure:

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: www.NeuLog.com. Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
 - a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

Tablet, smart phone with Bluetooth connectivity

Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

Procedure:

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

Operation with NeuLog Graphic Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

Procedure:

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

Off-line experiments

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

Materials needed:

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

Materials needed to configure your offline experiment:

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

Procedure:

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

Operation with NeuLog Digital Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

Procedure:

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

Note: If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

NeuLog Photo Gate Logger Sensor Guide NUL-209

NeuLog photo gate logger sensor NUL-209

The NeuLog photo gate sensor can be used for any science experiment which involves taking accurate velocity and/or acceleration measurements especially in the field of Physics.

The sensor comes pre-calibrated so you can start experimentation right out of the box using this guide.

Among hundreds of possible experimental subjects that can be studied with the NUL-209 sensor are: velocity/acceleration/distance relationships, impulse, collisions, conservation of energy, force, momentum, and many more.

The photo gate sensor uses an infrared beam which detects if an object is passing through it. Therefore, the sensor actively reads out as "open" or "blocked". While collecting data however, there are several parameters to truly customize your experiment:

- Velocity with a single gate
- Velocity with two gates
- Acceleration with a single gate
- Acceleration with two gates
- Change in time between two gates
- Velocities with timing card.

For the "two gates" options you will need two separate photo gate modules.

Experiment customization is explained after the following quick start guide.

Notes:

The photo gate sensor cannot work together with any other NeuLog sensor.

The photo gate sensor cannot work in the off-line mode since it is mandatory for it to be connected to the computer for its operation options. The off-line mode is disabled when the application finds that the photo gate sensor is connected.

The photo gate sensor cannot work with the Graphic display module, VIEW-101.


Experiment customization

The NeuLog photo gate sensor offers several different experimental customization options. To access them, click on the "Experiment Setup" icon after your photo gate sensor has been detected.


Timing cards:

Making timing cards is a unique option to further customize any experiment with the NeuLog photo gate sensor. To make timing cards we recommend cutting your desired shapes out of black plastic for the most accurate readings though even plain paper works just fine.


Timing cards examples:

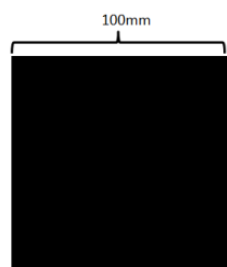For free fall experiments:


Single 'flag' timing card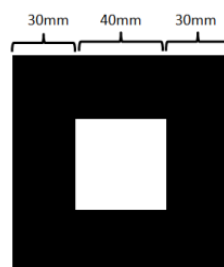


Double 'flag' timing card

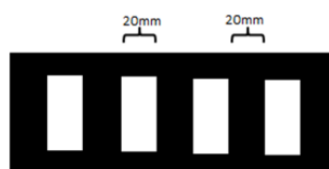Multi 'flag' timing card

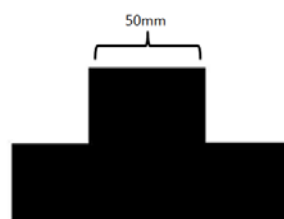For moving cart experiments:

Single 'flag' timing card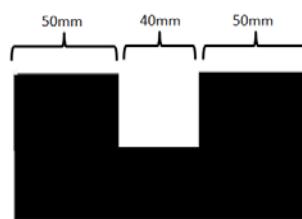

Double 'flag' timing card

Velocity with a single gate:

Velocity with a single gate is the very basic velocity measurement using only one photo gate sensor and any object with a known width.

1. Click the **Run Experiment** icon in the NeuLog application and select "Velocity with single gate".
2. The figure shows the basic idea of what your experimental setup should look like. An icon showing a blue photo gate sensor with a white beam and a small orange card passing through it.
3. Measure and input the object width (in millimeters) into the text field under the figure labeled "X (mm)".
4. Click on the **Record** icon.
5. For each time that the card passes through the photo gate, you will receive the time it took for it to pass and its velocity.

Acceleration with a single gate:

Acceleration with a single gate requires an object (or timing card) with two flags (example shown above) with known widths.

1. Click the **Run Experiment** icon in the NeuLog application and select "Acceleration with single gate".
2. The following figure shows the basic idea of what your experimental setup and timing card/object should look like.

3. Measure and input the width for both flags.
4. Click on the **Record** icon.
5. For each time that the card passes through the photo gate, you will receive its acceleration.

Acceleration with two gates:

Acceleration with two photo gates requires two NeuLog photo gate sensors and an object or timing card with a known width.

1. Click the **Run Experiment** icon in the NeuLog application and select "Acceleration with two gates".
2. The following figure shows the basic idea of what your experimental setup should look like.

3. Measure and input the width of your object (in millimeters) into the text field under the figure labeled "X (mm)".
4. Assign each of your photo gate sensors to both "Sensor A" and "Sensor B" position.
5. Click on the **Record** icon.
6. For each time that the card passes through the two photo gates, you will receive its acceleration.

Velocity with two gates:

The velocity with two gates option is ideal for experiments which examine collisions between two objects of known width and mass.

1. Click the **Run Experiment** icon in the NeuLog application and select "Velocity with two gates".
2. The following figure shows the basic idea of what your experimental setup should look like.

3. Measure and input the width and mass for both objects passing through the photo gates.
4. Assign each of your photo gate sensors to both "Sensor A" and "Sensor B" position.
5. Be sure to space the photo gates far enough apart to allow for a collision between the two objects before they rebound back through the photo gates.
6. Click on the **Record** icon.
7. For each time the cards pass through the photo gates, you will receive the velocity and momentum of each one of them.

Change in time between two gates:

Change in time between two photo gates requires two NeuLog photo gate sensors and an object or timing card with a known width.

1. Click the **Experiment Setup** icon in the NeuLog application and select "Delta between two gates".
2. The following figure shows the basic idea of what your experimental setup should look like.

3. Measure and input the width of your object (in millimeters) into the text field under the figure labeled "X (mm)".
4. Assign each of your photo gate sensors to both the "Sensor A" and "Sensor B" position.
5. Click on the **Record** icon.
6. For each time that the card passes through the two photo gates, you will receive the time difference between the first photo gate blocked and the second photo gate blocked.

Velocities with timing card:

Velocities with timing card experiments require one NeuLog photo gate sensor and an object or timing card (a multi 'flag' is recommended). A known width is not required for the time measurements, it is however required for velocity calculations.

1. Click the **Experiment Setup** icon in the NeuLog application and select "Velocities with timing card".
2. The following figure shows the basic idea of what your experimental setup should look like.

3. Choose the expected experiment. A longer duration will result in a smaller sample rate.
4. Click on the **Record** icon.
5. You will get the time difference in which every black flag passed through the photo gate. Velocities can be calculated by dividing the flag width by the time difference.

Included with sensor

- NeuLog General Guide.
- Infrared transmitter and receptor housed in the photo gate unit.
- Threaded rod for easy connection to retort stands or other lab equipment.

Sensor specifications	
ADC resolution	Digital
Accuracy	100 μ s
Resolution	100 μ s
Max sample rate (S/sec)	10,000

Sensor's features

- Fully digital data.
- Rugged plastic ergonomic case.
- Internally housed infrared transmitter and detector.
- Photo gate handle with two threaded holes to customize your experimental setup.
- Pushbutton switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).
- Pre-calibrated sensing equipment

Note: NeuLog products are intended for educational use.

Videos and experiment examples

- Videos, literature and other probes can be found at www.neulog.com.
- In order to access the photo gate sensor's page, choose "Products" on the main menu and then "Photo gate logger sensor".
- In order to find the access the photo gate sensor's experiments, choose "Example Labs":

Velocity of Free Falling Objects (P-35)

Horizontal Launched Projectile Motion (P-49)

Technical background

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (for example: °C, °F, Lux, %, ppm).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

The NeuLog photo gate sensor uses an infrared beam to determine when an object has passed through the photo gate. In every point where there is a break in the infrared beam, a 1 is registered in the program.

The NeuLog sensor firmware calculates velocities and accelerations based on the record of 1's and 0's as well as the experimental inputs. The computer program gets processed data from the sensor.

Maintenance and storage

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the photo gate sensor's body and probe.
- After use, gently wipe away any foreign material from the photo gate sensor.
- Store in a box at room temperature out of direct sunlight.

Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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