User Guide NUL-218

English

Español

NeuLog logger sensors and modules

NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

Note: NeuLog products are intended for educational use.

NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at www.NeuLog.com. Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

Note: There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to www.NeuLog.com before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

Quick start procedure

PC or Mac Computer – Wired Connection

Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at www.NeuLog.com as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

Procedure:

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

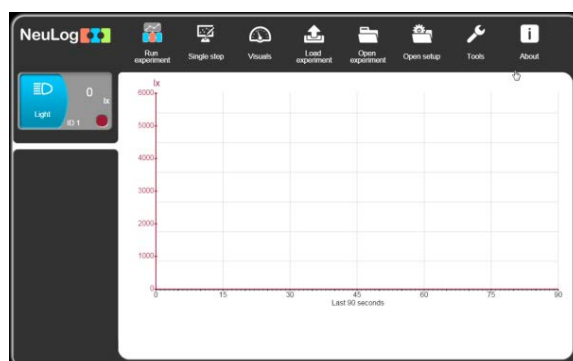6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- 'Run Experiment' with the connected sensors.
- 'Single Step' experiment.
- 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- 'Load Experiment' from the sensors' flash memory.
- 'Open Experiment' from a saved file.
- 'Tools' for changing the software settings or sensor ID.

Tablet, smart phone or any device with WIFI connectivity

Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

Procedure:

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: www.NeuLog.com. Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
 - a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

Tablet, smart phone with Bluetooth connectivity

Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

Procedure:

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

Operation with NeuLog Graphic Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

Procedure:

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

Off-line experiments

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

Materials needed:

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

Materials needed to configure your offline experiment:

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

Procedure:

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

Operation with NeuLog Digital Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

Procedure:

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

Note: If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

NeuLog Electrocardiogram Logger Sensor Guide NUL-218

NeuLog electrocardiogram logger sensor NUL-218

The NeuLog electrocardiogram sensor can be used for any science experiment that utilizes heartbeat data such as in the fields of: Biology, Physiology, Human medicine, Psychology, etc.

An electrocardiogram is a very commonly used medical device which measures and graphs the heart's electrical activity. Every single heartbeat is a very complex event which involves many anatomical locations undergoing changes in their electrical charges. A depolarizing cascade event originates in the Sinoatrial (SA) Node, often referred to as the natural pacemaker, and travels throughout the rest of the heart.

The sensor comes pre-calibrated so you can start experimentation right out of the box using this guide.

Among hundreds of possible experimental subjects that can be studied with the NUL-218 sensor are: electrical impulses, heart rate monitoring, human physiology, muscle physiology, heart rate changes to stimuli, and many more.

The ECG sensor's measurement unit is:


Arbitrary analog unit (Arb): An arbitrary unit to demonstrate waves, frequencies, and periods. What you see on the Y-axis is signal intensity, while the X-axis shows time.

Electrocardiograms

The heart undergoes a very complex process of electrical activity; this is the basis for how electrocardiograms work. As the heart's muscles contract, the cells are electrically depolarized causing a cascade of depolarization to flow across the heart – contracting the heart's muscle fibers.

During a heartbeat a complex series of events happen which begin with the cells resting potential at roughly -90 mV (millivolts). The heart's Sinoatrial Node begins the cascade by depolarizing – creating an electrical voltage. The electrical signal travels to the right atrium and almost immediately to the left atrium. Because the atria and the ventricles are insulated from each other by a special type of cell which acts as an electrical insulator, the signal is picked up in another cluster of cells called the Atrioventricular (AV) Node located in the right atria. The signal is carried by a special type of fibers – the Bundle of His into the ventricles. Purkinje fibers rapidly spread the electrical depolarization cascade throughout the ventricles. The entire event takes some time to complete. The graph is a visual demonstration of the heart's relative charge at any given point during heartbeats.

The electrical voltage is measured on the skin using conducting leads and it is converted into a graph that is very standard and reproducible.

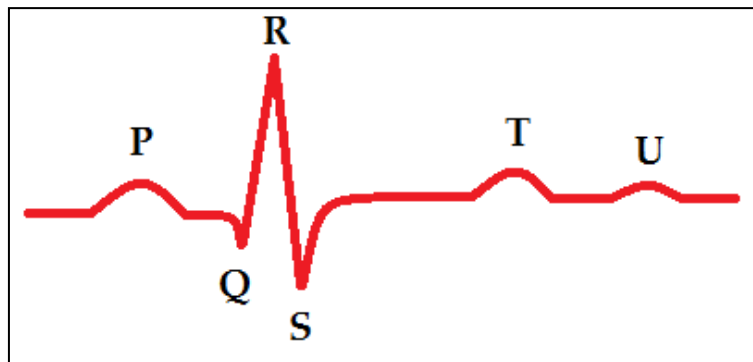- P – Contraction of atria
- Q – Initial depolarization of the interventricular septum
- R – Early depolarization of the ventricles
- S – Late depolarization of the ventricles
- T – Repolarization of the ventricles
- U – Often hidden by the T wave – this wave has an unknown origin

The graph in the previous page represents one complete heart beat from depolarization of the Sinoatrial Node to repolarization of the ventricles.

The following graph is from a NeuLog ECG experiment.

Connecting to the ECG sensor

The NeuLog electrocardiogram sensor uses three reusable clip leads for easy connection to the arms and legs. The heart's electrical fluctuations can be measured as a voltage signal appears between the left and the right arms.

This is a noisy signal and very difficult to measure. Instead, it is much better to use a reference electrode (usually connected to the left leg) and to measure the voltages of the left arm and the right arm in reference to this electrode and to amplify the difference between them. This is how it is done in NUL-218.

Leads:

Black – Reference
Red – Left
Yellow – Right

The NUL-218 is completely isolated from electrical power and signal in order to prevent the effect of these on the test subject.

How to connect:

1. Connect the NeuLog ECG sensor to a smart device using one of the following guides (PC or Mac/Tablet/Viewer).
2. Wet the red clip and place it on the subject's left wrist with the metal connector on the underside (palm side) of the arm.
3. Wet the yellow clip and place it on the subject's right wrist with the metal connector on the underside (palm side) of the arm.
4. Wet the black clip and place it on the subject's left leg or forearm or elbow.
5. Connect the three alligator clips (red, yellow, black) directly to their corresponding arm clip's (red, yellow, black) metal piece that forms a loop.
6. You are now connected.

Note: There are other configurations possible, this is just one. If the R wave shows a downward slope instead of upward, your connection's polarization is reversed though the absolute results are still the same.

Included with sensor

- NeuLog General Guide.
- NeuLog ECG probe with three alligator clips for easy connection to clamp leads.
- Three ECG clamp leads (red, yellow, black).

Sensor specifications	
Range and operation modes	0 to 4092 Arbitrary units
ADC resolution	12 bit
Resolution	1
Max sample rate (S/sec)	100

Experiment Duration: 1 second to 31 days.

Sensor's features

- Fully digital data.
- Rugged plastic ergonomic case.
- Pushbutton switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).
- ECG probe attached to the sensor's body by a durable rubber-coated wire with three alligator clip ends.
- Pre-calibrated sensing equipment.

Note: NeuLog products are intended for educational use.

Videos and experiment examples

- Videos, literature and other probes can be found at www.neulog.com.
- In order to access the electrocardiogram sensor's page, choose "Products" on the main menu and then "Electrocardiogram logger sensor".
- In order to access the ECG sensor's experiments, choose "Example Labs":

Electrocardiogram Measurement (B-40)

Technical background

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (for example: °C, °F, Lux, %, ppm).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

The ECG sensor detects minor changes in electrical voltage which originate in the heart and spread out across the body over a very short period of time. Cells have a resting potential of about 0.9 mV. A heart beat originates from an electrical impulse in the Sinoatrial Node, often called the natural pacemaker. The electrical potential cascade spreads across the heart. Using three leads (left, right and reference) the heart's electrical voltages (which originate in the heart) are measured.

Although the NeuLog ECG sensor is not aimed for medical measurements, it is built as an ECG real device.

One important feature is complete electrical isolation from the USB and the computer to which it is connected.

The power supply is isolated and transferred by transformers.

The signals are converted and transmitted by optical couplers.

Maintenance and storage

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the ECG sensor's body.
- After use, gently wipe away any foreign material from the ECG sensor.
- Store in a box at room temperature out of direct sunlight.

Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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