# User Guide NUL-219

English

Español

# NeuLog logger sensors and modules

## NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

**Note:** NeuLog products are intended for educational use.

## NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

**Note:** There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to [www.NeuLog.com](http://www.NeuLog.com) before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

## Quick start procedure

### PC or Mac Computer – Wired Connection

#### Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at [www.NeuLog.com](http://www.NeuLog.com) as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

## Procedure:

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

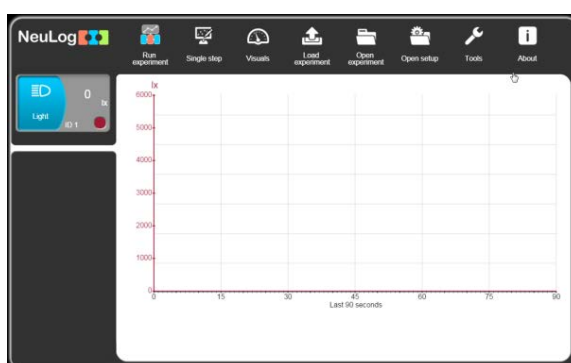6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- 'Run Experiment' with the connected sensors.
- 'Single Step' experiment.
- 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- 'Load Experiment' from the sensors' flash memory.
- 'Open Experiment' from a saved file.
- 'Tools' for changing the software settings or sensor ID.

### Tablet, smart phone or any device with WIFI connectivity

#### Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

**Procedure:**

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
  - a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

### Tablet, smart phone with Bluetooth connectivity

#### Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

**Procedure:**

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

### Operation with NeuLog Graphic Display Module (viewer)

#### Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

**Procedure:**

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

**Off-line experiments**

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

**Materials needed:**

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

**Materials needed to configure your offline experiment:**

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

**Procedure:**

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

## Operation with NeuLog Digital Display Module (viewer)

### Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

**Procedure:**

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

**Note:** If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

# NeuLog Colorimeter Logger Sensor Guide NUL-219

## NeuLog colorimeter logger sensor NUL-219

The colorimeter sensor can be used for any science experiment where liquid samples of different colors and concentrations are utilized such as in the fields of Chemistry, Biology, Pharmacology, Biochemistry, Environmental Science, Physics, or even just to study the behavior of light waves.

Colorimeters are used to determine absorbance or transmittance of given wavelengths (generally red, green, and blue) in a liquid solution. Absorbance and transmittance are directly related to the concentration of known solutes in the solution which is calculated using the Beer-Lambert law. One parameter can be calculated from the other.

The sensor comes pre-calibrated so you can start experimentation right out of the box using this guide.

Hundreds of possible experiments that can be performed with the NUL-219 sensor are: Analyzing chemical reactions, studying Beer's Law, water quality testing, monitoring of ecological systems, study of photosynthesis, studying light waves. The colorimeter could take the place of a spectrophotometer in many experiments.

The NeuLog colorimeter sensor's measurement units are:

- % Transmittance: The amount of light that passes through a sample.
- Absorbance: The amount of light which is absorbed by a sample.

## Cuvette usage and handling

It is important to maintain a standard for cuvette usage as their quality will directly impact the results of your experiments. If the cuvette has fingerprints, smudges, or cracks the sensor will detect them and give off an improper colorimeter reading. Smudges left from handling can easily be wiped off with tissue paper.

### Proper usage:

1. Carefully remove the cuvettes from their packaging.
2. Load your liquid sample into the cuvette before placing it into the sensor to prevent spilling and damaging the sensor.
3. Plug the cap into the cuvette and gently invert a few times. Shaking will create air bubbles which will be detected by the sensor and give an incorrect reading.
4. Before placing the cuvette in the sensor; hold it as far towards the top as you can and wipe off the bottom with a tissue. Do not touch the bottom of the cuvette again as this will leave a residue which could affect the results of the experiment.

5. If you have a cuvette with four clear sides, place it in the colorimeter sensor.

If you have a cuvette with two opaque sides:

- a. Determine which sides of the cuvette are clear and which sides have an opaque finish.
  - b. The clear opposing sides must be positioned facing the front and back of the sensor (they won't fit the other way) while the glossy sides face either side.
  - c. If the cuvette does not fit with a gentle push, rotate  $90^\circ$  in either direction and try again.
6. After experimentation, empty your samples, gently rinse out the cuvettes with distilled water and let them air-dry.

## Selecting a wavelength

Selecting the correct wavelength is important to get the best data and can easily be determined using one of two different methods. There is no incorrect wavelength, just some which are more informative than others. Lab experiments often come with a predetermined wavelength to use. If your experiment does not, follow one of these methods for best results:

- Knowing that the color we see solutions is based on the color reflected from them (and not absorbed): Choose a color you know will be absorbed. For example, if the solution is red try the blue LED option.
- Once the sensor is connected and the NeuLog software is open (follow one of the guides below for more instruction) click the “Module setup” button in the module box and try the different LED color options to determine which has the highest absorbance (it will appear in real time on the sensor’s module box). Select the color which has the highest absorbance reading.

### Included with sensor

- NeuLog General Guide.
- Three plastic cuvettes with three lids.

Sensors specifications		
	Transmittance	Absorbance
Range and operation modes	0-100% red, green, blue, and orange	0-4 red, green, blue, and orange
ADC resolution	14 bit	
Resolution	0.02%T	0.01 abs.
Max sample rate (S/sec)	100	

**Experiment Duration:** 1 second to 31 days.

### Sensor's features

- Fully digital data.
- Rugged plastic ergonomic case.
- Pushbutton switch for Start/Stop experiments in off line mode.
- Red, green, blue and orange LEDs for four different wavelength tests.
- Testing in both transmittance and absorbance
- Internal photodiode for detecting light.

**Note:** NeuLog products are intended for educational use.

## Videos and experiment examples

- Videos, literature and other probes can be found at [www.neulog.com](http://www.neulog.com).
- In order to access the colorimeter sensor's page, choose "Products" on the main menu and then "Colorimeter logger sensor".
- In order to access the colorimeter sensor's experiments, choose "Example Labs":

Beer-Lambert law (C-28)

Color Absorption (C-31)

## Technical background

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (for example: °C, °F, Lux, %, ppm).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

Inside the sensor's plastic body there are four LED lights (red, green, blue and orange) to test various wavelengths.

Light is transmitted from four different colored LEDs inside the sensor's body, one color at a time, passed through the cuvette containing a solution and then detected at the other end to determine the absorbance or transmittance of light at a specific wavelength.

The sensor works off of Beer's Law which is explained in the section "Beer's Law" below.

The four wavelengths emitted are:

- Red: 634-662 nm
- Blue: 452-476 nm
- Green: 500-530 nm
- Orange: 595-625 nm

### Beer's Law

Beer's Law is an extremely useful concept for solution concentration testing in the fields of Chemistry and Biology.

Standard curves are created by using solutions of known concentrations which are analyzed in a colorimeter to determine the absorbance and then plot a graph. The absorption for unknown samples can easily be tested and plotted on the standard curve to determine their concentration.

The law states that there is a logarithmic dependence between the transmission  $T$  of light through a substance, the product of the absorption coefficient of the substance  $\alpha$ , and the distance the light travels through the material  $l$ .

For liquids, these relations are usually written as:

$$T = \frac{I}{I_0} = 10^{-\alpha l}$$

Where:  $I_0$  and  $I$  are the intensity of the incident light and the transmitted light, respectively.

The transmission (or transmittance) is expressed in terms of absorbance which, for liquids, is defined as:

$$A = -\log_{10} \left( \frac{I}{I_0} \right) = -\log_{10}(T)$$

This implies that the absorbance becomes linear with the concentration (or number density of absorbers) according to:

$$A = \alpha l$$

## Maintenance and storage

- Never submerge the NeuLog plastic body in any liquid.
- Never place anything in the colorimeter sensor's opening on the faceplate except the approved cuvettes.
- Do not allow liquid into the NeuLog plastic body.
- After using the probe wipe off all excess material, liquid or residue from the colorimeter probe.
- Store in a box at room temperature out of direct sunlight.
- The cuvette handling and storage guide can be found at the beginning of this document.

## Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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