User Guide NUL-240

English

Español

NeuLog logger sensors and modules

NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

Note: NeuLog products are intended for educational use.

NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at www.NeuLog.com. Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

Note: There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to www.NeuLog.com before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

Quick start procedure

PC or Mac Computer – Wired Connection

Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at www.NeuLog.com as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

Procedure:

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

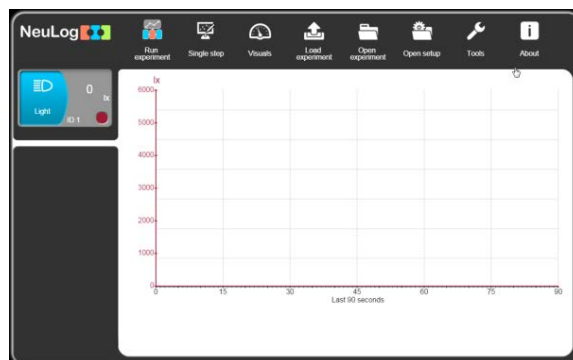6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- a. 'Run Experiment' with the connected sensors.
- b. 'Single Step' experiment.
- c. 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- d. 'Load Experiment' from the sensors' flash memory.
- e. 'Open Experiment' from a saved file.
- f. 'Tools' for changing the software settings or sensor ID.

Tablet, smart phone or any device with WIFI connectivity

Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

Procedure:

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: www.NeuLog.com. Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
 - a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

Tablet, smart phone with Bluetooth connectivity

Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

Procedure:

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
 - b. **'Single Step'** experiment.
 - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
 - d. **'Load Experiment'** from the sensors' flash memory.
 - e. **'Open Experiment'** from a saved file.
 - f. **'Tools'** for changing the program settings or sensor ID.

Operation with NeuLog Graphic Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

Procedure:

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

Off-line experiments

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

Materials needed:

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

Materials needed to configure your offline experiment:

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

Procedure:

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

Operation with NeuLog Digital Display Module (viewer)

Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

Procedure:

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

Note: If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

NeuLog Ammonium Logger Sensor Guide NUL-240

NeuLog ammonium logger sensor NUL-240

The NeuLog ammonium ion sensor can be used for any science experiment which utilizes ammonium concentration readings such as in the fields of Environmental Science, Ecology, Chemistry, Biology, etc.

Among hundreds of possible experimental subjects that can be studied with the NUL-240 sensor are: chemical reactions, water quality, environmental health, small animal behavior, ecological studies, and many more.

The ammonium sensor's measurement units are:

- Milligrams per liter (mg/L): SI measurement unit for density of a solid dissolved in a liquid.
- Parts per million (ppm): The total amount of a specific molecule (ammonium for example) per million molecules in a sample.

Ammonium sensor usage

The ammonium sensor's electrode uses a PermaFil (non-refillable), ion-selective electrode. The reference chamber is filled with gel and sealed; therefore, no reference filling solution is required.

Usage guide:

| | |
|-----------------------------|--------------------------------------------------|
| Concentration range | 1M to 5×10^{-6} M 0.01 to 18,000 ppm |
| pH range | 4 to 10 pH |
| Temperature range | 0 to 50°C |
| Minimum sample size | 3mL in a 50 mL beaker |
| Reproducibility | ±4% |
| Electrode resistance | 1 to 4 megaohms |

Equipment required:

- Wash bottle with distilled or deionized water.
- Several clean beakers.
- 1 mL, 10 mL and 50 mL pipettes (a 50 ml graduated cylinder can also be used).

Technical notes:

- All standard solutions should be prepared fresh.
- The standard solutions should be at the same temperature as the sample.
- Use ISA in all solutions (also recommended for the measured sample).

Note: ISA is added to all solutions to ensure the samples and standard solutions have the same ionic strength; the ratio is 1:50.

Electrode preparation

1. Remove the protective cap encasing the ammonium probe's electrode.

CAUTION: Do not touch the PVC membrane at the end of the probe with your fingers.

2. Rinse the electrode with deionized water and blot dry. Be sure not to rub as this will harm the sensor.
3. Soak the electrode in deionized water for 10 minutes, then in a diluted ammonium standard solution (0.0001 M or 1 ppm) for two hours until ready for use.
4. For preparing the 1 ppm standard solution, mix one drop of the 1000 ppm standard solution with 20 drops of ISA and add deionized water up to a total of 40 ml solution (assuming one drop equals 0.04 ml). It is also possible to use a micropipette if available.

Offsetting the sensor

Be sure to offset the ammonium sensor once daily during periods of regular usage.

1. Prepare the electrode as described in the "Electrode preparation" guide above. If you have already prepared the electrode in the last three days and stored it as described in the "Electrode storage" section then you do not have to repeat it before offsetting the sensor.
2. Prepare the offsetting solution: Mix 1 ml of the ISA solution with 0.5 ml of the 1000 ppm standard solution. Add deionized water up to a total of 50 ml. You will get a 10 ppm solution.
3. Rinse the electrode with deionized water, blot dry and place in the beaker.
4. After a stable reading has been reached, press the "Start/Stop" button located on the sensor's faceplate for 4 seconds.
5. Rinse the electrode once more with deionized water and blot dry.
6. The ammonium sensor is now ready for use.

Electrode storage

Short term (overnight or weekend):

Rinse the electrode thoroughly with deionized water and place the tip in a diluted standard solution (around 10 ppm) between measurements.

Long term:

Rinse the electrode thoroughly with deionized water and blot dry. Attach the hard plastic protective electrode cap again to protect the sensing element. Follow procedures in the "Electrode preparation" section when using the electrode again.

| |
|----------------------|
| Included with sensor |
|----------------------|

- NeuLog General Guide.
- NeuLog ammonium (NH_4^+) sensor electrode.
- Ammonium (NH_4^+) Ionic Strength Adjuster (ISA).
- Ammonium (NH_4^+) 1000 ppm standard.

| Sensor specifications | | |
|---------------------------|------------------------------------------------------------------------------------|---------------------------------------------------------------------------------|
| | mg/L | ppm |
| Range and operation modes | 0.02 to 18,000 | 0.02 to 18,000 |
| ADC resolution | 15 bit | |
| Resolution | 0.04 at 0.02 to 100 mg/L 0.4 at 100 to 1,000 mg/L 40 at 1,000 to 18,000 mg/L | 0.04 at 0.02 to 100 ppm 0.4 at 100 to 1,000 ppm 40 at 1,000 to 18,000 ppm |
| Max sample rate (S/sec) | 100 | |

Experiment Duration: 1 second to 31 days.

Sensor's features

- Fully digital data.
- Rugged plastic ergonomic case.
- PermaFil ion selective electrode.
- Ammonium probe connected to the sensor's body by a flexible rubber-coated wire.
- Protective hard-plastic storage cap.
- Pushbutton switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).

Note: NeuLog products are intended for educational use.

Videos and experiment examples

- Videos, literature and other probes can be found at www.neulog.com.
- In order to access the ammonium sensor's page, choose "Products" on the main menu and then "Ammonium Logger Sensor".
- In order to access the ammonium sensor's experiments, choose "Example Labs":

Ammonium and Nitrate in an Aquarium (B-45)

Technical background

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (for example: °C, °F, Lux, %, ppm).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

The ammonium sensor uses a permanently filled ion selective electrode to accurately measure the amount of ammonium ion in a solution.

The Ionic Strength Adjustor (ISA) is used to standardize the strength of the ions in various solutions to give the most accurate readings.

Maintenance and storage

Before use, please review the "Ammonium sensor usage", "Electrode preparation", "Electrode storage", and "Offsetting the sensor" sections of this document above to ensure the proper storage and longevity of the NeuLog ammonium ion sensor.

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the ammonium sensor's body.
- After use, gently wipe away any foreign material from the ammonium sensor.
- Store in a box at room temperature out of direct sunlight.

Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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