# User Guide PANDA-1

English

Español

# NeuLog logger sensors and modules

## NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

**Note:** NeuLog products are intended for educational use.

## NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

**Note:** There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to [www.NeuLog.com](http://www.NeuLog.com) before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

## Quick start procedure

### PC or Mac Computer – Wired Connection

#### Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at [www.NeuLog.com](http://www.NeuLog.com) as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

**Procedure:**

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

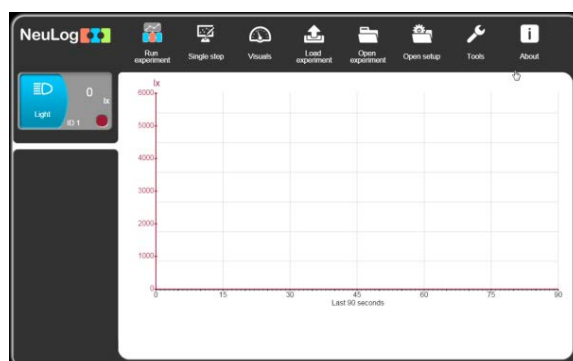6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- 'Run Experiment' with the connected sensors.
- 'Single Step' experiment.
- 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- 'Load Experiment' from the sensors' flash memory.
- 'Open Experiment' from a saved file.
- 'Tools' for changing the software settings or sensor ID.

### Tablet, smart phone or any device with WIFI connectivity

#### Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

**Procedure:**

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
  - a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

### Tablet, smart phone with Bluetooth connectivity

#### Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

**Procedure:**

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

### Operation with NeuLog Graphic Display Module (viewer)

#### Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

**Procedure:**

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

**Off-line experiments**

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

**Materials needed:**

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

**Materials needed to configure your offline experiment:**

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

**Procedure:**

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

## Operation with NeuLog Digital Display Module (viewer)

### Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

**Procedure:**

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

**Note:** If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

# NeuLog Panda Multi-Sensor Guide

## NeuLog Panda multi-sensor PANDA-1

The NeuLog Panda multi-sensor can be used for any science experiment which utilizes Room Temperature, Light, Barometric pressure, Altitude, Sound, Relative Humidity, Dew-Point, Magnetic Field and G-Acceleration readings.

Panda is used for environmental data collection and other experiments in Natural Science, Biology, Physics and Chemistry.

The multi-sensor comes pre-calibrated so you can start experimentation right out of the box using this guide.

Hundreds of possible experimental subjects can be performed with the Panda multi-sensor. Examples are: Light, relative humidity and temperature at different environments, Qualitative and quantitative description of nature, Sound intensity levels, Sound and distance, Rate of heating and cooling, Barometric pressure and altitude measurements at different heights, Barometric pressure and weather, Magnetism – positive and negative poles, Compass experiments, Acceleration in an elevator or car, Acceleration and harmonic motion, Dew point and dew formation.

The internal sensors' measurement units are:

### Temperature sensor:

- **Celsius (°C):** The SI (International System of Units) unit of temperature.
- **Fahrenheit (°F):** The temperature measurement unit of the English System commonly used in the United States.

### Light sensor:

The measurement unit for all four data collection ranges (low, medium, high) is the lux. Lux (lx, or lux): The SI unit of light intensity

### Barometer sensor:

The barometer sensor can collect data using four different units of measurement:

- **Kilopascal (kPa):** The SI unit of pressure
- **Atmospheres (atm):** A pressure unit equal to one Earth atmosphere at sea level (101.325 kPa)
- **Inches of mercury (in Hg):** A non-SI unit of atmospheric pressure
- **Millimeters of mercury (mm Hg):** A non-SI unit of atmospheric pressure

### Altitude sensor:

The altitude sensor's units are Meters (M), the SI base unit of length.

### Sound sensor:

The sound sensor's measurement unit is:

**Decibel (dB):** A measurement unit that shows intensity (loudness of sound). This is a logarithmic unit.

### Relative humidity sensor:

Data is presented as a percentage which is the ratio of the amount of water vapor in the air versus the maximum amount that could be allowed at a given temperature.

### Dew point sensor:

The NeuLog dew point temperature is presented in the following units of measurement:

- **Celsius:** The SI (International System of Units) unit of temperature.
- **Fahrenheit:** The temperature measurement unit of the English System commonly used in the United States.

### Magnetic field sensor:

The magnetic field sensor's measurement units are:

- **Micro Tesla ( $\mu\text{T}$ ):** The SI unit of magnetic flux density (magnetic field).
- Angle degrees ( $^{\circ}$ ).

### G-Acceleration sensor:

The acceleration sensor's measurement unit is:

**Meters per second squared ( $\text{m/s}^2$ ):** The SI unit of acceleration or change in velocity over time.

The Panda enables to zero offset the G-acceleration value in order to turn it to acceleration sensor.

Included with sensor
----------------------

- NeuLog General Guide.
- USB to mini USB cable.
- 9 Internal sensors.
- Internal rechargeable battery that can be charged by the cable (battery level is shown on screen).

Sensor specifications		
Range and operation modes	ADC resolution	Resolution
<b>Temperature sensor</b>		
-25 to 60 °C	12 bit	0.1 °C
-13 to 140 °F	12 bit	0.2 °F
<b>Light sensor</b>		
1,000 lx	16 bit	1 lx
6,000 lx	16 bit	1 lx
60,000 lx	16 bit	1 lx
240,000 lx	18 bit	6 lx
<b>Barometer sensor</b>		
66 to 110 kPa	24 bit	0.1 kPa
0.65 to 1.08 atm	24 bit	0.01 atm
19.49 to 32.48 in Hg	24 bit	0.01 in Hg
495 to 825 mm Hg	24 bit	0.1 mm Hg
<b>Altitude sensor</b>		
-722 to 3000 m	24 bit	0.9 m
<b>Sound sensor</b>		
40 to 110 dB	12 bit	0.1 dB
<b>Humidity sensor</b>		
0 to 100 %	16 bit	0.1 %


Sensor specifications		
Range and operation modes	ADC resolution	Resolution
<b>Dew point sensor</b>		
-114 to 109°C	12 bit	0.1 °C
-182 to 228°F	12 bit	0.2 °F
<b>Magnetic field sensor</b>		
X Axis -200 to 200 μT	16 bit	0.1 μT
Y Axis -200 to 200 μT	16 bit	0.1 μT
Z Axis -200 to 200 μT	16 bit	0.1 μT
0° to 360°	16 bit	1°
<b>Acceleration sensor</b>		
X Axis -20 to 20 m/s <sup>2</sup>	16 bit	0.01 m/s <sup>2</sup>
Y Axis -20 to 20 m/s <sup>2</sup>	16 bit	0.01 m/s <sup>2</sup>
Z Axis -20 to 20 m/s <sup>2</sup>	16 bit	0.01 m/s <sup>2</sup>

NeuLog Panda can work with external sensors (up to 5 at a time) as well (these are not included in the Panda package).

Part No.	Sensor
NUL-201	Voltage
NUL-202	Current
NUL-203	Temperature
NUL-204	Light
NUL-205	Oxygen
NUL-206	pH sensor
NUL-208	Heart rate & pulse
NUL-209	Photo gate
NUL-210	Pressure
NUL-211	Force
NUL-212	Sound
NUL-213	Motion
NUL-215	Conductivity
NUL-218	Electrocardiogram
NUL-219	Colorimeter
NUL-220	CO <sub>2</sub>
NUL-222	Blood pressure
NUL-230	UVB
NUL-232	UVA
NUL-243	GPS position

## Panda's features

- Fully digital data.
- Rugged plastic case.
- Pushbutton switch that powers the multi-sensor on and off.
- Sensor icons that show the internal sensors.

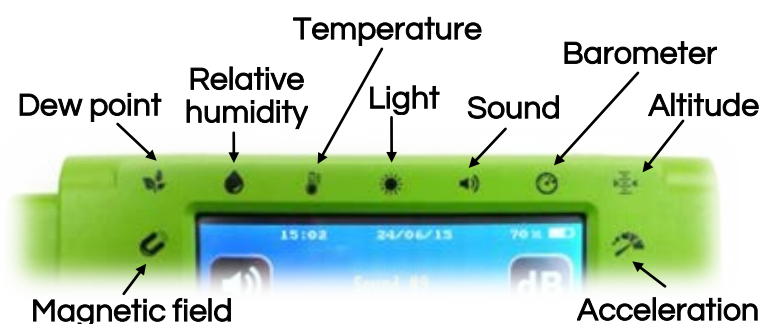- Sensors are pre-calibrated.
- 3.2" (320X240 pixels) Color display touch screen that turns off after 2 minutes without use (4 minutes until full device turns off).
- Has the ability to save 20 experiments in the internal memory (no matter what the experiment parameters are).

**Note:** NeuLog products are intended for educational use.

## Technical background

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This is also true for the Panda Multi-sensor. This technology allows the multi-sensor to collect and then store the digital data in the correct scientific units (for example: °C, °F, Lux, %).

The multi-sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using provided firmware update.

The viewer's screen uses resistive touchscreen technology which reacts to a pressure being applied to the screen – opposed to the more common capacitive screen which requires an electrical conductor to operate. This allows for students to operate the screen even while wearing gloves.

### Temperature:

Inside the Panda device is a thermistor which is a variable resistor that varies significantly with temperature. The resistance increases nonlinearly with a decreasing temperature. During calibration, the sensor's internal controller creates the conversion formula from resistance to temperature.

## Light:

Inside the hard plastic casing, lies a photodiode which reacts with photons to release free electrons (photoelectrons). The amount of light striking the sensor is directly proportional to the voltage generated by the photoelectrons released. The sensor measures the general voltage released and thus calculates the light intensity.

If the light readout is very low, try changing the sensor's mode to a higher sensitivity.

## Barometer:

The barometer sensor uses the piezoresistive effect. The transducer is made of two metal foils separated by silicon; when pressure is applied on the transducer its resistance changes. One side of it is at complete vacuum which enables the measurement of the absolute pressure on its other side.

## Altitude:

The altitude reading is based on the barometric pressure which is translated to meters.

## Sound:

The sound sensor uses an internal microphone with a series of circuits, filters, and amplifiers to best isolate a specific sound source. Sound waves enter through the hole at the top of the sensor's plastic body so you should point it directly towards the sound source for best readings.

**Relative humidity:**

The relative humidity sensor's casing houses a capacitive polymer which reacts with water vapor present in the air. The reaction with the polymer produces a voltage which is relative to the concentration of water vapor in the air.

**Dew point:**

Dew point is calculated using both the environmental temperature and relative humidity values.

**Magnetic field:**

The magnetic field sensor utilizes the Hall Effect to generate a voltage relative to the magnetic field. The Hall Effect is a natural phenomenon where a magnetic field exerts a force on charge carriers if an electrical current flows through the magnetic field. The charge differential is measured and can easily be converted into the magnetic field strength because of the direct relationship.

**G-Acceleration:**

The G-acceleration IC sensor houses three small thin silicon accelerometers. Each accelerometer is oriented along a different axis (X, Y, or Z) connected to a weight and based on piezoresistive technology. When an accelerometer experiences force along its axis, the silicon changes its resistance proportionally to the force it is experiencing.

These resistor changes are converted into voltage by a Wheatstone bridge circuit and read by the IC internal controller. The IC converts the force into acceleration and transmits, upon request, the 3D accelerations in a digital form.

## Making measurements with NeuLog Panda

Panda is operated by independent internal firmware for performing experiments and data collection.

The Panda firmware is specially designed for elementary school level; it is unique, super friendly and very intuitive.

In **normal mode**, Panda displays the value measured of one of its internal or external sensors in one display option:


In **experiment record** mode, Panda displays the sampled values of up to two sensors at a time. It records, in its flash memory, the measurements of all its internal sensors (and also external if connected while the measurement is performed). These values can be displayed at any time.

## Application

Panda's application (for PC, MAC or tablets) has the same features as the internal Panda software along with the following ones:

1. Full view option (displays all 4 options together):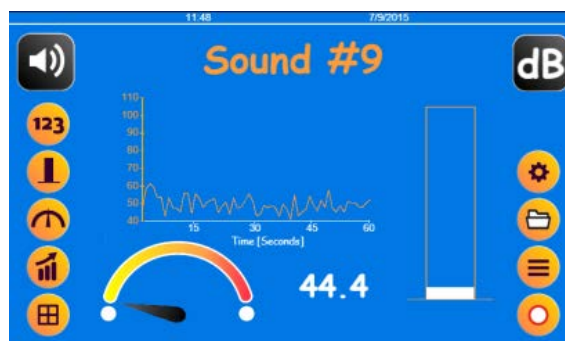


2. Saving experiments as CSV files after recording.
3. Opening CSV files.
4. After recording an experiment, a cursor can be added to focus a certain point.
5. Tools: Sync date and time (synchronizes the data and time from the computer to the top of the screen).

## Acceleration sensor orientation

The NeuLog acceleration sensor (an accelerometer) measures static acceleration due to gravity and dynamic acceleration due to the sensor's motion. By measuring static acceleration, one can find out the angle the sensor is tilted with respect to the Earth. By measuring dynamic acceleration, one can find out how the velocity of the sensor changes with time.

When the sensor is not moving and is facing upwards, the reading is  $9.8 \text{ m/s}^2$  (because it includes static acceleration). When the sensor is at free fall and is facing upwards, the reading is  $0 \text{ m/s}^2$ . If we are interested only in the vertical axis (Z axis) and in the dynamic acceleration, we can zero it by pressing the zero button. This zero offset is canceled when turning the Panda off and on.

The Panda acceleration sensor has the ability to measure acceleration in three different axes (X, Y and Z). The figure below demonstrates the axes.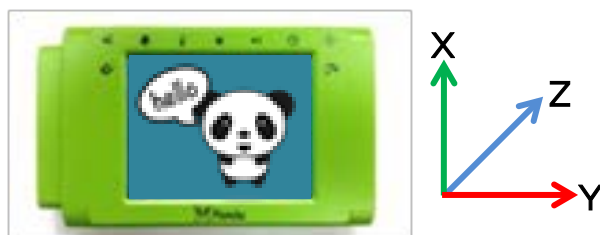


## Magnetic sensor orientation

When pointing the positive side to east or west, the value should be close to  $0 \mu\text{T}$ .

When pointing the positive side to the north, the value should be close to  $30 \mu\text{T}$ .

When pointing the positive side to the south, the value should be close to  $-30 \mu\text{T}$ .

The Panda magnetic field sensor has the ability to measure magnetic field in three different axes (X, Y and Z). The figure below demonstrates the axes.

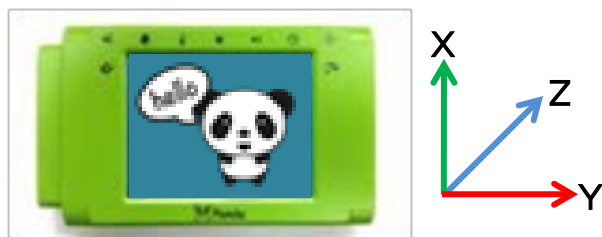It is important to work in an environment with minimal metals, since they can influence the sensor's value.

## Using Panda with WiFi

Materials needed:

- NeuLog Panda multi-sensor
- WIFI-201 or WIFI-202 WiFi Module

Your Panda Multi-sensor needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing (tablet, smartphone, or computer). Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

**Procedure (follow these steps in this exact order):**

1. The WiFi module should be off at this point (not connected to anything).
2. Take you panda and press on the settings icon (on the right side of the Panda screen).
3. Press on WiFi mode.
4. Connect the Panda multi-sensor directly to the WiFi module (no wires required).
5. Although not required, we recommend plugging the Panda to an outlet using a USB to mini USB charger (such as a typical cell phone charger).

6. The WiFi module indicator lights will flash; take no action until the LED to the far left turns blue; this can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for the example).
8. Give your device about 20 seconds to connect to the WiFi module.
9. Once the device is connected, go to your browser and type the website [panda201.com](http://panda201.com) into the URL bar, then wait for 30-60 seconds.
10. The Panda multi-sensor will give a live reading and you can begin experimentation.

## Maintenance and storage

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the Panda multi-sensor body.
- After use, gently wipe away any foreign material from the sensor.
- Store in a box at room temperature out of direct sunlight.

## Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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