# User Guide WIFI-202

English

Español

# NeuLog logger sensors and modules

## NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

**Note:** NeuLog products are intended for educational use.

## NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

**Note:** There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to [www.NeuLog.com](http://www.NeuLog.com) before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

## Quick start procedure

### PC or Mac Computer – Wired Connection

#### Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at [www.NeuLog.com](http://www.NeuLog.com) as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

**Procedure:**

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

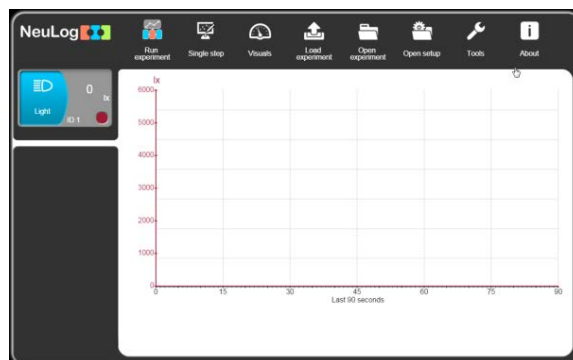6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- 'Run Experiment' with the connected sensors.
- 'Single Step' experiment.
- 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- 'Load Experiment' from the sensors' flash memory.
- 'Open Experiment' from a saved file.
- 'Tools' for changing the software settings or sensor ID.

### Tablet, smart phone or any device with WIFI connectivity

#### Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

**Procedure:**

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
  - a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

### Tablet, smart phone with Bluetooth connectivity

#### Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

**Procedure:**

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

### Operation with NeuLog Graphic Display Module (viewer)

#### Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

**Procedure:**

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

**Off-line experiments**

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

**Materials needed:**

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

**Materials needed to configure your offline experiment:**

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

**Procedure:**

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

## Operation with NeuLog Digital Display Module (viewer)

### Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

**Procedure:**

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

**Note:** If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

# NeuLog WIFI Module Guide WIFI-202

## NeuLog WIFI module WIFI-202

The WIFI module is without a doubt one of the most advanced and innovative elements that NeuLog has to offer. Data collection and analysis with NeuLog sensors can be performed through any device which has wireless capabilities such as iPads, Android tablets, smart phones, Windows, Mac, and Linux based computers. The WIFI module can seamlessly transform group projects by connecting up to 5 devices onto the wireless network at a single time – promoting interaction and allowing each member to have their own dataset.

There is no downloaded application or software installation as they are built directly into the WIFI module which transmits its own closed wireless network. This means that you do not need your own wireless network – though if you have one, a simple setting can be changed on the WIFI module to allow surfing of the web at the same time!

All current popular browsing programs can be used with the WIFI module to allow you to use whichever browser you and your students feel most comfortable with.

## Using the NeuLog WIFI module

The NeuLog WIFI module is a very advanced piece of equipment; having capability to transmit its own wireless network as well record data from up to 5 NeuLog sensors at a time.

The NeuLog WIFI module can be powered by the NeuLog battery module or by a USB power source.

It also can be used as a USB module when it is connected to a PC or Mac and under local NeuLog software.

You'll notice that on the front of the WIFI-202 module there is a LED indicator that includes four color LEDs. Underneath the indicator, there is a legend that explains what each color represents:

- **Blue (access point)** – When the blue LED is the only light on, the WIFI module is in Access Point mode (described below) and is transmitting a closed wireless network which you can connect devices to.
- **Green (client mode)** – When the green LED is the only light on, the WIFI module is in Client Mode (described below) and allows your wireless devices to control sensors through the website [www.wifi201.com](http://www.wifi201.com) as well as browse the internet.
- **Red (transmission)** – The red LED will blink when there is active communication between any connected sensors and the computer or smart device.
- **Orange (USB module)** – The orange LED will remain lit when the WIFI module is working in USB mode (detailed below). USB mode can be turned on and off by pressing the button on the front of the WIFI module three times.

## Materials required

- Battery module or USB to micro USB cord (as a power source)
- Any NeuLog sensor
- NeuLog WIFI module
- Any device with wireless capabilities.

## Control and view modes

When multiple users are connected to the WIFI module's network; one user at a time can enter "control mode" the others users will be defaulted to "view mode".

### Control mode:

Entering control mode allows that user to manipulate the experimental parameters and begin data collection using the web application – this prevents multiple users from all trying to change the experiment at once.

To leave the control mode, simply click the "Tools" icon and then the "Release master" button.

### View mode:

In view mode; users have almost full access to the web application – only lacking the ability to manipulate data and start/stop experiments. Users in view mode have the full ability to view data in real time.

**Important:**

In the event that anyone used the WIFI module as the controller and forgot to disconnect from it, the WIFI module will not allow any other user to connect to the WIFI module as a controller until it is restarted.

In order to allow another controller to access the WIFI module, press and hold the push button located on the front panel of the WIFI module for 3 seconds until all 3 lights turn on – this will refresh the module resetting it back to having no users in "Control Mode".

**Access point mode use**

Access point mode is the default operating mode for the the NeuLog WIFI module – it transmits its own closed wireless network which allows up to 5 devices to connect to it at a time. No prior wireless networks are required, making this option both very portable and unique. Because the network is closed; you will not be able to browse the internet, if you wish to browse the internet while using the WIFI module please locate the "Client Mode" procedural guide in this manual. Connect up to 5 NeuLog sensors directly to the left side of the WIFI-202 module (no wires required).

1. Connect either a BAT-200 or a USB to micro USB cable connected to a power source, to the WIFI-202 module to supply power.
2. The WIFI-202 indicator lights will flash; take no action until the LED to the far left turns blue.
3. Take your tablet or smart phone and go to the WiFi settings and select the NeuLog network which matches the WIFI module ID found on the back of the WIFI-202 device (NeuLog1334 for example).
4. Once the device is connected go to your browser and type in the website [wifi201.com](http://wifi201.com) into the URL bar, then wait for 30-60 seconds.

**Note:** Only one device can be in control mode at a time, the remaining devices must be in "View mode" but switching between them is both possible and very easy. The "Control mode" allows the user to change experimental parameters and both start and stop data collection. View mode allows the user to view the data being recorded.

5. The browser will then load a new screen and begin to auto detect the sensors, this can take a minute.
6. Once the sensors are found you will see an icon on the left side of the screen for any detected sensors. The icon will display data in real time.
7. To set the sensors' settings click on each sensor's icon.

8. To change the experimental settings click the "Run experiment" icon (experiment duration for example).
9. To run an experiment and collect data click "Record".

### Client mode use

If you have a wireless network and wish to browse the internet while using the NeuLog WIFI module; you can use client mode to connect the WIFI module to your local internet. By default the WIFI module starts in "Access point mode" but can easily be changed to "Client mode".

**Note:** Before following this procedural guide you must configure your WIFI module using the "Access Point Mode" guide above. Access point mode is most commonly used.

1. Configure your WIFI module as per the "Access point mode" guide in this document.
2. Using your connected device, locate and click the "Tools" icon on the top of the NeuLog web application window in your browser.
3. Click on "WIFI options".
4. Input your local internet connection's name and password in their respective text fields and click on "Connect". This step attempts to connect the NeuLog WIFI module directly to your local wireless network allowing you to browse the internet.

**Note:**

During this time the green light on the WIFI module will flash. In the event that the NeuLog site could not be accessed the blue light will turn back on and the WIFI module will return to "Access point mode".

5. If connection to the NeuLog site is successful; you will be disconnected from the NeuLog closed wireless network (i.e. NeuLog1334) and the green light on the WIFI module will remain on.
6. At this point you must connect to your local wireless connection (the information you put into the "Client Mode" boxes).
7. On the new screen you get, click the Connect button to reconnect the browser to the WIFI-202 module.
8. You will now be in client mode with the ability to browse the internet using your smart device.

When connecting additional wireless devices (up to 5 total) simply type the address "wifi201.com" into your device's internet browser bar. A prompt will ask for the WIFI-202 ID, this is the digit code located on the back of the WIFI module. Click "Connect" and now you will be able to access the NeuLog software broadcast from the WIFI module.

The next time the WIFI module is turned on it will again attempt to enter client mode by connecting to the same wireless connection. If the connection fails; the WIFI module will restart back in access point mode.

## Using the WIFI-202 in USB mode

USB mode allows you to use your WIFI module as if it were a NeuLog USB-200. This requires a computer which has the NeuLog software (available as a free download on [www.NeuLog.com](http://www.NeuLog.com)).

By connecting the WIFI module to your sensors and computer via the included USB to micro USB cable, you are able to stream data directly from your sensors to the NeuLog browser based application.

### To enter USB Mode:

1. Download and install the NeuLog browser based application onto the computer which you plan on using. The application can be downloaded for free at [www.NeuLog.com](http://www.NeuLog.com) as well as a full application user guide.
2. Choose "Downloads" on the main menu and then choose "Software and Application".
3. Connect the WIFI module to the computer with the NeuLog application using the included USB to micro USB cable, you'll see the LED lights blink for a moment while the module powers on.
4. Connect your NeuLog sensors to the side of the WIFI module, the red LED on each sensor should blink to acknowledge they are working properly.

5. Press the button located on the front panel of the WIFI module three times, the yellow LED labeled 'U' should turn on.
6. Open the NeuLog application.
7. Your connected sensors will automatically be detected.
8. Once your sensors are detected, you can begin experimenting in USB mode!

**Note:** To switch back to Access Point mode from USB mode, press the button on the WIFI module faceplate three times again.

### Troubleshooting walkthrough

Due to the technological complexity of the NeuLog WIFI-202 module and the ever updating software and hardware found on today's wireless devices, it is possible that you may run into some performance issues while the unit is communicating with your wireless device. Luckily, for these circumstances the NeuLog WIFI module can easily be cleared and updated!

If you are noticing that the performance of your NeuLog WIFI-202 module is acting sporadic please follow the guides below.

## Restarting the WIFI module:

Restarting the WIFI module will often clear up any connectivity issues you may be having and is an extremely easy process.

1. Connect the WIFI module to a power source – this can be either a charged BAT-200 or directly to your computer using the USB to micro USB cable.
2. The LED lights on the WIFI module will turn on to indicate that it is now being powered.

**Note:** You do NOT need to wait for the network to be transmitted at this point.

3. Once the WIFI module is powered, simply press and hold the button on the front panel of the WIFI module for roughly 4 seconds (until the green, red, and blue turn on).
4. When the blue LED is the only light on, the closed wireless network is being transmitted and you can connect devices as usual.
5. If the behavior of the WIFI module is still sporadic please continue to the **Updating the WIFI module's firmware** walkthrough below.

## Updating your WIFI module's firmware:

If restarting your NeuLog WIFI module didn't solve the WIFI module's performance issues, please update your WIFI module firmware as per the guide below.

Firmware updates for the NeuLog WIFI module are regularly released to the NeuLog website [www.NeuLog.com](http://www.NeuLog.com). Choose "Downloads" on the main menu and then choose "Upgrade firmware".

It is recommended to check back often for the latest releases.

**Note:** A full WIFI firmware update walkthrough is also available on the website listed above.

### Materials required

- A computer with wireless capability running a Windows operating system
- WIFI-202 module + USB to micro USB cable

1. Download the latest NeuLog WIFI firmware update from the NeuLog website.
2. Install the NeuLog WIFI firmware updater.
3. Power your WIFI module by connecting a charged BAT-200 or through connection to your computer using the included USB to micro USB cable.
4. After a few moments when only the blue LED light remains on – the wireless network (NeuLog1334 for example) is being transmitted.
5. Open your computers wireless connection settings and connect to the NeuLogXXXX wireless network.

**Note:** You may get a notification that you now have a "limited" connection – this is OK as it simply means you don't have access to the internet.

6. Open the WIFI firmware updater, titled "wifi201 update" from your desktop (the file you downloaded and installed in step 1).
7. A small window will open. Press Start. The updater will run automatically and will take about five minutes.

8. When the firmware update is complete (you will see the word "Status: Update finish OK" at the bottom of the window) you can close the window, and power cycle the WIFI module by unplugging it from your power source.
9. Your WIFI module is now up to date!
10. You can verify it by checking the version on the "About NeuLog" icon (after entering [wifi201.com](http://wifi201.com)).

If problems with your NeuLog WIFI-201 module continue to persist, please contact NeuLog for additional assistance at [info@neulog.com](mailto:info@neulog.com).

#### Included in package

- NeuLog General Guide.
- NeuLog WIFI-202 module.
- USB to micro USB cable.

#### Maintenance and storage

- Never submerge the NeuLog plastic body in any liquid.
- Store in a box at room temperature out of direct sunlight.

## Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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